

Technical Information

Indumax H CLS54

Inductive conductivity sensor with hygienic design for applications in the food, beverage and pharmaceutical industries and in biotechnology







Standard Number 74-03

With quality certificate

Application

The CLS54 conductivity sensor is specifically intended for use in hygienic applications in the food, beverage and pharmaceutical industries and in biotechnology. Thanks to its food-grade virgin PEEK body, its jointless and creviceless design and its hygienic certificates, it meets the extremely exacting demands of these industries. The CLS54 is especially suitable for the following applications:

- Phase separation of product/water and product/product mixtures in pipe systems
- Control of CIP (cleaning in place) processes in the return channel
- Concentration control in remaking of CIP cleaning agents
- Product monitoring in pipe systems, bottling plants, quality assurance
- Leakage monitoring
- in the following industries
- Dairies
- Breweries
- Beverages (water, juices, soft drinks)
- Pharmaceuticals and biotechnology

The sensor can be used with the transmitters Liquiline M CM42, Liquisys M CLM223/253 and Mycom S CLM153; it is part of the measuring system Smartec S CLD134 and compatible with the transmitter electronics of Smartec S CLD132.

Your benefits

- Unique hygienic design, thus no risk of recontamination
- With all hygienic certificates required in hygienic applications
- With all process connections commonly used in hygienic applications
- Durable thanks to encapsulated, jointless design
- Fast measurement with temperature response time t₉₀ under 26 s enables safe and efficient phase separations
- Insensitive to electrode soiling and polarization thanks to electrodeless, inductive measuring principle
- Suitable for application in hazardous areas



People for Process Automation


TI400C/07/en/07.10 71108219

Function and system design

Measuring principle

Inductive conductivity measurement

A generator (1) generates an alternating magnetic field in the primary coil (2) which induces a current in the medium (3). The strength of the induced current depends on the conductivity and thus the ion concentration of the medium. The current flow in the medium generates another magnetic field in the secondary coil (4). The resulting current induced in the coil is measured by the receiver (5) and processed to determine the conductivity.

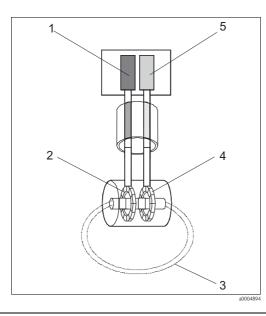

Inductive conductivity measurement

- 1 Generator
- 2 Primary coil
- 3 Current flow in the medium
- 4 Secondary coil
- 5 Receiver

Benefits of inductive conductivity measurement

- No electrodes, therefore no polarization
- Accurate measurement in media or solutions with a high soiling degree and a tendency to deposition
- Complete galvanic separation of measurement and medium

Important properties Indumax H CLS54

Hygiene

The injection-molded sensor is made of highly chemically, mechanically and thermally resistant PEEK. It does not have joints or crevices and is therefore hygienically safe. The exclusive use of virgin PEEK as material in contact with medium guarantees highest biological safety for the food, beverage and pharmaceutical areas. The sensor has been designed according to the effective guidelines of ASME BPE (The American Society of Mechanical Engineers – Bioprocessing Equipment).

Certified hygiene

The sensor has all necessary listings and approvals for hygienic areas, such as exclusive use of FDA-listed materials for surfaces in contact with medium and certification according to EHEDG guidelines (European Hygienic Engineering and Design Group) as well as 3-A. A certificate on the biological reactivity test according to USP part <87> and <88> class VI including lot tracing back to compounding is available as option.

Process connections

The sensor is available with all process connections commonly used in hygienic applications. The most common connections are available as standard. Further process connections can be ordered as special versions.

Process temperature, process pressure

The use of special components and materials makes the sensor suitable for continuous exposure to temperatures of +125 °C (275 °F). Short-time (max. 60 min.), it may be exposed to +150 °C (300 °F) for sterilization. The sensor is pressure-resistant up to 12 bar (174 psi) up to 90 °C (194 °F). At higher temperatures, its pressure resistance is always higher than the respective steam pressure. The sensor is suitable for underpressure applications.


Temperature measurement

The sensor has an integrated temperature sensor with a temperature response time $t_{90} < 26$ s. This enables economically efficient phase separations at varying and quickly changing process temperatures. The temperature sensor is embedded into the PEEK body eliminating the need for seals which ensures a long service life.

Measuring system

A complete measuring system comprises:

- a CLS54 inductive conductivity sensor
- a transmitter, e.g. Liquiline CM42

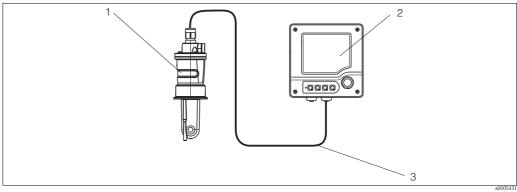

BN n.c.

Measuring system example

- Indumax H CLS54 1
- 2 3 Transmitter Liquiline CM42
- Measuring cable

Input

Measured values	Conductivity Temperature		
Cell constant k	$k = 6.3 \text{ cm}^{-1}$		
Measuring ranges	Conductivity Temperature	recommended range: 100 μ S/cm to -10 to +150 °C (+14 to +302 °F)	2000 mS/cm (uncompensated)
Temperature measurement	Pt 1000 (class A acc. to IEC 6	50751)	
Cable specification	Indumax H is supplied with a special measuring cable.	fixed cable. The connection to the transmitte	er can be extended using the CLK5
		Cable	
	v	NH Coax WH { Ferrule RD Ferrule BU Coax RD { Ferrule RD Ferrule BU Ferrule BU	Screen Inner connector Screen Inner connector
		GN GN KWH YE	

Fixed cable / CLK5 special measuring cable

Cable length: max. 55 m (180 ft) total length


Performance characteristics

Temperature response time	$t_{90} \le 26 \text{ s}$
Measured error	\pm (0.5 % of measured value + 10 $\mu S/cm)$ after calibration (plus inaccuracy of the conductivity of the calibration solution)

Installation

Installation instructions

The sensor has to be immersed completely into the media. Avoid bubbles in the area of the sensor.

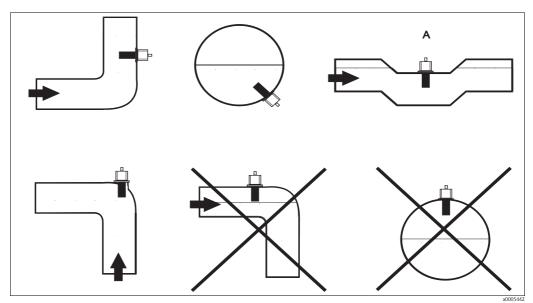

Installation positions of conductivity sensors

A Installation position is not usable for hygienic applications.

The media should flow through the flow opening of the sensor (see indicator arrow of the sensor body). The symmetrical measuring channel allows a flow in both directions.


In narrow installation conditions, the ion flow in the medium is affected by the pipe walls. This effect is compensated by the so-called installation factor. The installation factor can be entered in the transmitter or the cell constant can be corrected by multiplication with the installation factor to ensure correct measurement.

The value of the installation factor depends on the diameter and the conductivity of the pipe as well as the sensor's distance from the wall.

If the distance from the wall is sufficient (a > 15 mm, from DN 65), it is not necessary to consider the installation factor (f = 1.00).

If the distance from the wall is smaller, the installation factor increases in case of electrically insulating pipes (f > 1) and decreases in case of electrically conductive pipes (f < 1).


The installation factor can be measured using calibration solutions or it can be approximately determined from the diagram beside.

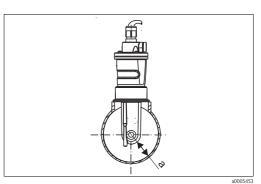



а

Sensor distance from the pipe wall

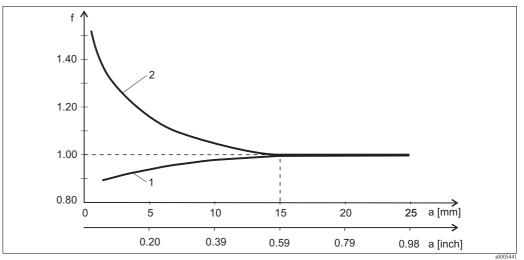

Dependance of installation factor f on wall distance a

1 Conductive pipe

2 Insulating pipe

Air set

To compensate residual coupling in the cable and between the two sensor coils, you must perform a zero calibration in air ("air set") before installing the sensor. For further information, refer to the Operating Instructions of your transmitter.

Environment

Ambient temperature	-20 to +60 °C (-4 to +140 °F)
Storage temperature	-25 to +80 °C (-13 to +176 °F)
Relative humidity	5 to 95 %
Ingress protection	IP 68 / NEMA 6P (1m (3.28 ft)water column, 50 °C (122 °F), 168 h)


	Process			
Process temperature	-10 to +125 °C (+14 to +257 °F)			
Sterilization	150 °C / 5 bar (302 °F / 72.5 psi) (max. 60 min.)			
Process pressure	12 bar (174 psi) up to 90 °C (194 °F) 8 bar (116 psi) at 125 °C (257 °F) 0 to 5 bar (0 to 72.5 psi) when in CNR-applicable areas (tested with 50 bar (725 psi)) underpressure down to 0.1 bar (1.45 psi) absolute			
Pressure / temperature load curve	$\begin{bmatrix} psi \\ 174 \\ 12 \\ 116 \\ 8 \\ 72.5 \\ 58 \\ 4 \\ -10 \\ 0 \\ 10 \\ 30 \\ 50 \\ 70 \\ 90 \\ 110 \\ 125 \\ 150 \\ 10 \\ 125 \\ 150 \\ 10 \\ 10 \\ 125 \\ 150 \\ 10 \\ 10 \\ 125 \\ 150 \\ 10 \\ 10 \\ 10 \\ 125 \\ 150 \\ 10 \\ 10 \\ 10 \\ 10 \\ 10 \\ 10 \\ $			

Pressure / temperature load curve

- Α
- short-time sterilization (max. 60 min) MAWP (maximum allowable working pressure) according to ASME-BPVC Sec. VIII, Div 1, UG101 for CRN В registration


Mechanical construction

Design, dimensions

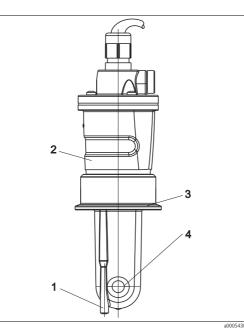

Indumax H CLS54

- 1 Temperature sensor
 - Housing
- 2 3 Process connection
- Flow opening 4

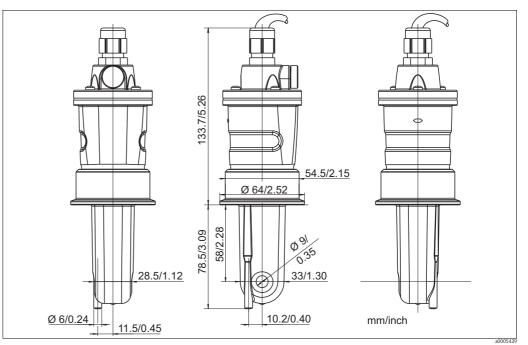

Dimensions CLS54 (long version)


Weight	Depending on version 0.3 to 0.5 k	Depending on version 0.3 to 0.5 kg (0.66 to 1.1 lb.) plus cable		
Surface roughness	$Ra \leq 0.8~\mu m$ (smooth, injection-mo	$Ra \leq 0.8~\mu\text{m}$ (smooth, injection-molded PEEK surface) for the surfaces in contact with medium		
Materials	In contact with medium Not in contact with medium	Virgin PEEK PPS-GF40 Stainless steel 1.4404 (AISI 316L) Screws: 1.4301 (AISI 304) PVDF (cable gland) FKM, EPDM (seal) TPE (cable)		

Process connections

- Dairy pipe fitting DIN 11851, DN 50 ^{a)}
- Aseptic fitting DIN 11864-1 form A, for pipe according to DIN 11850, DN 50
- Clamp ISO 2852 (also for TriClamp, DIN 32676), 2"
- SMS 2" b)
- Varivent N DN 40 125
- NEUMO BioControl D50 (DN 40.2")

Additional process connections are available on request.

- a) Dairy pipe fitting DIN 11851 is generally not considered hygienic. With the adapter SKS Siersma, it meets the 3-A standards requirements.
- b) Process connection is not considered hygienic according to the requirements of EHEDG.

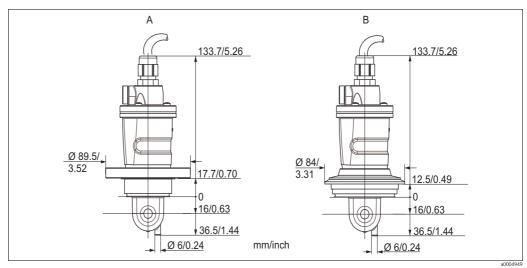


Process connections CLS54 (short version)

A NEUMO BioControl D50

for pipe connection: DN 40 (DIN 11866 series A, DIN 11850) DN 42.4 (DIN 11866 series B, DIN EN ISO 1127) 2" (DIN 11866 series C, ASME-BPE)

B Varivent N DN 40 to 125

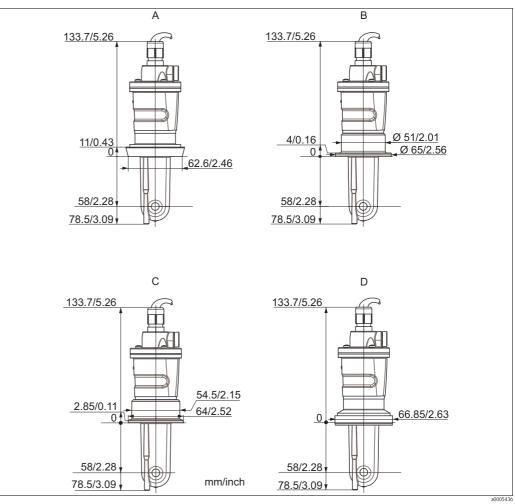

Process connections CLS54 (long version)

Dairy pipe fitting DIN 11851, DN 50 SMS 2" Α

В

- Clamp ISO 2852 , 2" С
- D Aseptic fitting DIN 11864-1 form A, for pipe according to DIN 11850, DN 50

Electrical connection

Fixed cable connection with cable gland

Chemical durability

Medium	Concentration	РЕЕК
Caustic soda NaOH	0 to 15 %	20 to 90 °C (68 to 194 °F)
Nitric acid HNO ₃	0 to 25 %	20 to 90 °C (68 to 194 °F)
Phosphoric acid H ₃ PO ₄	0 to 15 %	20 to 80 °C (68 to 176 °F)
Sulphuric acid H_2SO_4	0 to 30 %	20 °C (68 °F)
Peracetic acid H ₃ C-CO-OOH	0.2 %	20 °C (68 °F)

No responsibility is taken for the correctness of this information.

		All materials in contact with medium are listet at FDA.		
		EHEDG Certified for cleanability according to EHEDG document 2 Note! The cleanability of a sensor also depends on the way of installation. To install the sensor in a pipe system use		
		the appropriate and EHEDG certified flow assembly for the respective process connection. 3-A Certified according to 3-A Standard 74-03 ("3-A Sanitary Standards for Sensor and Sensor Fittings and Connections Used on Milk and Milk Products Equipment").		
		Biological reactivity (USP class VI) (optional) Certificate (Certificate of Compliance) on biological reactivity tests according to USP (United States Pharmacopeia) part <87> and part <88> class VI with traceability of the materials in contact with medium.		
Ex approvals		 ATEX II 1G Ex ia IIC T3/T4/T6 CSA IS/NI Cl. I Div. 1 & 2 GP A - D in connection with the transmitter Liquiline M CM42 		
Pressure approvals		Canadian pressure approval for pipes according to ASME B31.3		

Certificates and approvals

FDA

Ordering information

A	pproval	proval		
A	Non-E	Non-Ex area		
G	ATEX	II 1G Ex i	a IIC T3/T4/1	Γ6
0	CSA I	CSA IS/NI Cl. I Div. 1 & 2 GP A - D		
	Proc	ess conn	ection	
	MV5	Dairy pi	pe fitting DIN	11851, DN 50 ¹⁾
	AA5	Aseptic	fitting DIN 11	864-1 form A, pipe DIN 11850, DN 50
	CS1	Clamp I	SO 2852, 2" (1	ong)
	SMS			
	VA4	Varivent	t® N DN 40 to	125
	BC5	C5 NEUMO BioControl [®] D50		
		Supple	ementary ec	quipment
		0	Standard	
		2	Biological read	tivity tests according to USP <87>, <88> class VI
		3	CRN approval	(ASME B31.3) ³⁾
		4	CRN approval	$(ASME \ B31.3)^{3)}$ + biological reactivity tests according to USP <87>, <88> class V
			Cable lengt	h
			1 Fixed	cable 5 m (16.41 ft.)
			2 Fixed	cable 10 m (32.81 ft.)
			3 Fixed	cable 20 m (65.62 ft.)
			Tem	perature sensor
			2	Integrated temperature sensor Pt 1000
		1 1		complete order code

the 3-A standards requirements.

²⁾ Process connection is not considered hygienic according to the requirements of EHEDG.

³⁾ CRN approval only valid for process connections MV5, CS1 and VA4.

Hygienic certificates

Product structure Indumax H CLS54

	ACCOSULCS	
Measuring cables	 CLK5 measuring cable Extension cable for connecting CLS52/54 and transmitter via the VBM junction box, sold by the meter; order no. 50085473 	
	Junction box VBM For cable extension 10 terminals Cable entries: 2 x Pg 13.5 or 2 x NPT ½" Material: aluminum Ingress protection: IP 65 (≙ NEMA 4X) Order numbers: – cable entries Pg 13.5: 50003987 – cable entries NPT ½": 51500177	
Calibration solutions	 Precision solutions, traceable to SRM (standard reference material) by NIST, for qualified calibration of conductivity measurement systems according to ISO 9000, with temperature table CLY11-B 149.6 µS/cm (reference temperature 25 °C / 77 °F), 500 ml / 16.9 fl.oz Order no. 50081903 CLY11-C 1.406 mS/cm (reference temperature 25 °C / 77 °F), 500 ml / 16.9 fl.oz Order no. 50081904 CLY11-D 12.64 mS/cm (reference temperature 25 °C / 77 °F), 500 ml / 16.9 fl.oz Order no. 50081905 CLY11-E 107.0 mS/cm (reference temperature 25 °C / 77 °F), 500 ml / 16.9 fl.oz Order no. 50081905 	
Related products	 Indumax P CLS50 Highly resistant inductive conductivity sensor for standard, hazardous and high-temperature applications; for ordering information, see Technical Information TI182C/07/en Smartec S CLD134 Measuring system for conductivity and concentration measurement in the food industry; for ordering information, see Technical Information TI401C/07/en 	

Accessories

Documentation

Transmitters	 Mycom S CLM153, Technical Information TI234C/07/en Liquisys M CLM223/253, Technical Information TI193C/07/en Liquisys M CLM223F, Technical Information TI237C/07/en Liquiline M CM42, Technical Information TI381C/07/en 		
	 CLS54-G, Safety instructions for electrical equipment in hazardous areas, XA400C/07/a3 		
Calibration solutions	 Precision calibration solution CLY11, Technical Information TI162C/07/en 		

Instruments International

Endress+Hauser Instruments International AG Kaegenstrasse 2 4153 Reinach Switzerland

Tel.+41 61 715 81 00 Fax+41 61 715 25 00 www.endress.com info@ii.endress.com

71108219



People for Process Automation

